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The following pages were translated using Google Translate. For this reason, their correctness cannot be guaranteed. The German version of the operating instructions is legally binding.

Structural unit/working area:
Quantum optics

Activity: Working on laboratory fume hoods / digestors (standard laboratory fume hoods DIN EN 14 175)

DESIGNATION

fume hoods (standard laboratory fume hoods DIN EN 14 175)

DANGERS TO PEOPLE AND THE ENVIRONMENT

dangers to humans

- Danger due to escape of hazardous substances when the front lid is open, heavy construction, currents and turbulence in front of the exhaust, large thermal loads.
- Risk of explosion if large quantities of flammable gases, vapors, dust or aerosols are released.
- Danger of splashing of substances or of splinters and fragments being thrown out when the front lid is open.
- Danger of the front lid falling down if the cable breaks.

PROTECTIVE MEASURES AND RULES OF CONDUCT

Organizational protective measures

- Keep the front lid closed and work through horizontal slides or access openings if possible.
- Only place the material needed for experimentation into the fume hood when working.
- Do not place any bulky objects in the fume hood.
- Do not walk quickly past the trigger and do not lean into the trigger.
- Do not overload the fume cupboard with large releases of substances (gases, vapours); wherever possible and reasonable, collect and remove emissions at the point of release.
- Do not operate large thermal loads in the fume cupboard (do not operate several gas burners or several heating plates in one fume cupboard: otherwise use a smoke extractor (according to DIN 12924-2: 2007)).
- No storage of chemicals in the fume hood; this is intended for the safe conduct of experiments and is not a collection point for chemicals that are no longer needed.
- Compliance with workplace exposure limits (see **TRGS 900**) is the responsibility of the respective user. If necessary, the user must carry out/commission orientation measurements.

Personal protective measures

- As with all work in the laboratory, protective goggles must be worn when working in the fume hood.

BEHAVIOR IN THE EVENT OF DISTURBANCES

disorders

- If the exhaust air alarm sounds (horn, red light), do not continue working under any circumstances: Danger of hazardous substance release, danger of explosion – report to control room (Tel. No. **22225**).
- If you hear unusual noises, the front lid is stiff or tilted, stop working - report to the control room (Tel. No. **22225**).

ACCIDENTS - FIRST AID - EMERGENCY CALL 112


- Make an emergency call.
- Initiate first aid measures.
- Remove unconscious persons to fresh air while protecting yourself. Keep warm and calm. If breathing stops, give artificial respiration.
- Trained first aiders: see first aid notice
- First aid services must be recorded in the first aid book.

TESTING, MAINTENANCE AND REPAIR


maintenance and repair

- Inspections, maintenance and repairs are organized by Dept. V-2 (please register deductions there)
- Repairs may only be carried out by authorized persons.
- Properly tested fume cupboards carry a sticker (1), fume cupboards that operate with reduced performance are marked with (2). Here, the user must check and document the extent to which the functional limitation affects the safety of the employees.

(1)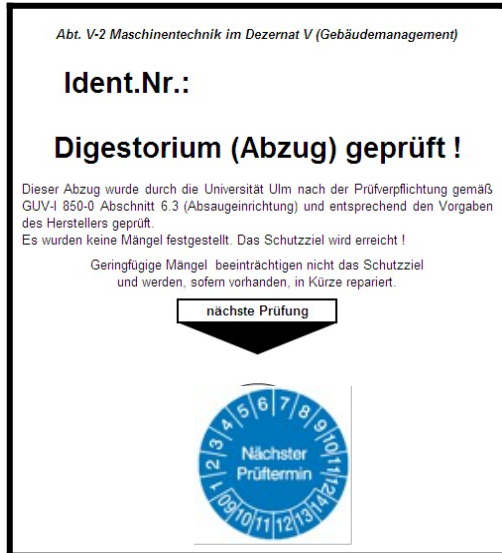

(2)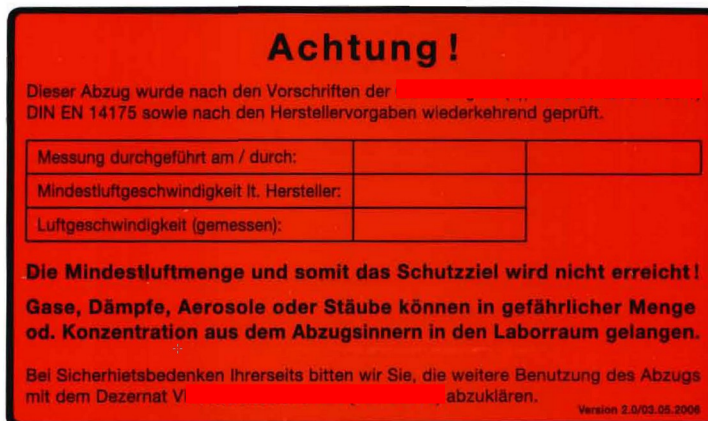

CONSEQUENCES OF NON-COMPLIANCE

injuries

- Improper handling can lead to injuries

property damage

- Misconduct can lead to contamination of the laboratory by hazardous substances.

Legal consequences

- Operating instructions are binding and represent written occupational safety instructions for employees.
- Failure to comply may result in legal consequences.
- Failure to follow this instruction is a violation of the given instructions and will be punished accordingly.

Creation date: 10.12.2024 – Creator: Michael Ferner