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The following pages were translated using Google Translate. For this reason, their correctness cannot be guaranteed. The German version of the operating instructions is legally binding.

# operating instructions

organizational unit: Quantum Optics

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## DESIGNATION

### Handling compressed gas cylinders in safety cabinets.

Basically, TRGS 510 applies to the proper storage of compressed gas cylinders.

## DANGERS TO PEOPLE AND THE ENVIRONMENT

- Formation of ignitable and/or explosive vapor-air mixtures in the usable space of the cabinet due to leaky gas bottles, fittings or pipes and insufficient extraction.
- Accumulation of suffocating gases in the usable space of the cabinet due to leaky gas bottles, fittings or pipes and insufficient extraction.

## PROTECTIVE MEASURES AND RULES OF CONDUCT

- The procurement, storage, use and necessary protective measures for pressure vessels must be discussed in advance with the safety officer.
- Only the designated gas cylinders may be connected and operated in the safety cabinets; care must be taken to ensure that the correct pressure reducer is used for the type of gas.
- The manufacturer's operating instructions must be observed.
- The installation and connection of gas cylinders may only be carried out by authorized and trained personnel.
- Only trained persons may open the safety cabinet.
- Safety cabinets may only be operated without restrictions if they are earthed, connected to the exhaust air and inspected by a qualified person.
- When opening the safety cabinet and when handling items in the cabinet, appropriate protective clothing, including safety goggles and a lab coat, must be worn.
- The cabinet surface must be kept clear at all times to avoid accidentally covering the air supply opening.
- The doors of safety cabinets must not be held open by wedges or objects placed in front of them, nor must they be locked when open.
- Dust deposits and other dirt in the cabinet must be removed regularly, otherwise the exhaust air performance will be reduced.

### Bottle change:

- To change the bottle, close the valve of the empty bottle, screw on the protective cap and remove the bottle from the cabinet.
- Place the full bottle in the cupboard and secure it with the holder.
- Connect the fittings in the bottle cabinet to the bottle, observing the fitting manufacturer's operating instructions.
- After opening the cylinder valve, check the connections for leaks using leak detection spray. Repair any leaks found.
- Close the cabinet door.
- When changing bottles, appropriate protective clothing, including safety shoes, must be worn.

## BEHAVIOR IN THE EVENT OF DISTURBANCES/IRREGULARITIES

- Repairs to the safety cabinet and fittings may only be carried out by qualified persons.
- If the cabinet ventilation fails, gas extraction from cylinders containing flammable and toxic/very toxic gases must be stopped immediately.
- Close the bottle valves on each bottle.
- Inform superiors.
- At a temperature of 50 °C the automatic closing mechanism is triggered. Do not reach into the cabinet again!
- After a fire, the safety cabinet may only be opened by the fire department.

FIRE DEPARTMENT / RESCUE CONTROL CENTER: 112  
Control room in case of technical faults: 22225

## ACCIDENTS - FIRST AID - EMERGENCY CALL 112

Where possible, treat injuries immediately using standard first aid measures and inform first responders.

Everyone should find out where the first aid kit and fire extinguisher are located at their workplace.

Report all accidents (even seemingly minor ones) to your supervisor and document them in the first aid report pad

Observe posted fire safety regulations and rescue plans; read them regularly so that no unnecessary time is lost in an emergency.

FIRE DEPARTMENT / RESCUE CONTROL CENTER: 112

COMPANY DOCTOR: 66190

## INSPECTION AND MAINTENANCE

### Monthly functional test by the user and before/when changing gas bottles:

- Checking the proper functioning of the doors including hinges, locking systems, door closers, door holding systems
- Visible defects e.g. on the cables or the ramp etc.

### Annual maintenance by a qualified technician:

- Safety cabinets for compressed gas cylinders must be checked regularly by a qualified person. Particular attention must be paid to the locking devices for doors and connections, the seals and the air exchange
- For inspection date, see inspection sticker
- Carry out maintenance work only after consultation with the responsible laboratory management

## CONSEQUENCES OF NON-COMPLIANCE

### health consequences

- Contact with gases can cause injury
- Gas leakage can lead to explosions
- Gas leakage can lead to suffocation

### property damage

- e.g. by explosions

### Legal consequences

- Operating instructions are binding and represent written occupational safety instructions for employees.
- Failure to comply may result in legal consequences.

Failure to follow this instruction is a violation of the given instructions and will be punished accordingly.