BEZEICHNUNG

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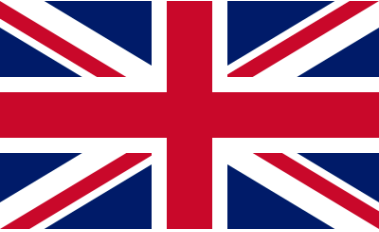
The following pages were translated using Google Translate. For this reason, their correctness cannot be guaranteed. The German version of the operating instructions is legally binding.

Structural unit/working area:
Quantum optics

Activity: Working with table centrifuges

DESIGNATION

Handling table centrifuges

DANGERS FOR HUMANS AND THE ENVIRONMENT

Dangers to humans

- Danger of rotor breakage if the rotor is not properly maintained!
- Release of dangerous substances and/or formation of dangerous aerosols if liquids leak or if rotors or centrifuge buckets break!
- Risk of fire and explosion when handling flammable substances!
- Danger of flying parts, e.g. if the rotor breaks!
- Hand and finger injury from contact with moving rotor

PROTECTIVE MEASURES AND RULES OF CONDUCT

Technical protective measures

- Do not remove or tamper with protective devices.
- Choose the working height of the centrifuge so that the rotor chamber is visible!
- Detection of vessel breakage and leakage is only possible with centrifuges with a transparent lid.
- The centrifuge must not be opened or interfered with during operation!
- Regularly clean the rotors according to the manufacturer's instructions! The cleaning must be documented, stating the agents used!
- Only use rotors and centrifuge cups/tubes that are approved for use in this centrifuge (see operating instructions)!
- Permanently mark the direction of rotation.
- The centrifuge must be placed securely on a suitable flat surface and a free space of at least 30 cm must be maintained around the centrifuge (GUV-R 500 and manufacturer's recommendations).

Organisational protective measures

- Only trained employees may be entrusted with the operation.
- Before operation, the rotor, centrifuge cups/tubes and their closures must be visually inspected for corrosion and integrity. Sealing rings must also be checked for integrity and proper seating!
- Centrifuges with defects must not be operated!
- The permissible load and speed must not be exceeded (see operating instructions). If it is absolutely necessary to exceed the permissible load, the volume or speed must be reduced in accordance with the operating instructions!
- Observe operating instructions/safety data sheets for the hazardous substances used.
- Rotors must be loaded evenly to prevent unacceptable imbalance and to minimize the risk of fluid leakage!
- After starting the centrifuge, employees should remain at the device until the set rotation speed is reached. During the waiting phase, it can be determined whether there is an imbalance or whether unclear noises on the device indicate a fault. In such cases, the centrifuge can be switched off immediately with an EMERGENCY STOP.

Personal protective measures

- When cleaning the system, it is recommended to wear protective goggles and, if necessary, gloves (only when the machine is at a standstill).

RESPONSE TO MALFUNCTIONS

Disturbances

- In the event of a malfunction, the device must be shut down immediately and - secured against being switched on again accidentally. The supervisor must be informed immediately.
- Malfunctions may only be eliminated to the extent instructed in accordance with the workplace-specific instructions.

Troubleshooting, repair and maintenance

- Before carrying out any maintenance or repair, always unplug the appliance.
- Portable electrical equipment must be inspected by a qualified expert at least once a year. The test results must be recorded in a test log.
- A test must be carried out in disassembled condition every three years. Exceptions to these test intervals can be made if the centrifuge is not explosive, flammable or explosive and the kinetic energy is less than 10,000 Nm or the rated power is less than 500 W (manufacturer's information).
- Repairs may only be carried out by authorized and trained personnel .

MAINTENANCE AND REPAIR

Maintenance and repair

- Maintenance and repair work may only be carried out when the machine is at a standstill.
- The device must be secured against unauthorized switching on.
- Repairs, maintenance work and inspections may only be carried out by authorized and qualified persons.

WHAT TO DO IN THE EVENT OF ACCIDENTS - FIRST AID - EMERGENCY CALL 112

- Make an emergency call.
- Initiate first aid measures.
- Trained first aiders: see first aid notice
- First aid services must be recorded in the first aid book.
- Bangs and explosions – even if no personal injury occurs – must be reported to the responsible occupational safety officer and the UKBW.

CONSEQUENCES OF NON-COMPLIANCE

Injuries

- Flying splinters/parts can cause injuries (especially to the eyes)
- Released hazardous substances can cause health damage

Property damage

- Misconduct may result in device failure.

Legal consequences

- Operating instructions are binding and represent written occupational safety - instructions for employees.
- Failure to comply may result in legal consequences.
- Failure to follow this instruction is a violation of the given instructions and will be punished accordingly.