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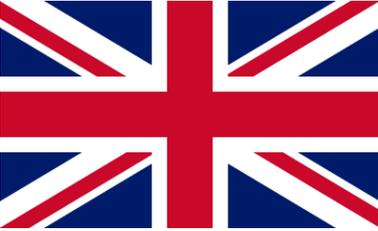
The following pages were translated using Google Translate. For this reason, their correctness cannot be guaranteed. The German version of the operating instructions is legally binding.

Structural unit/working area:
Quantum optics

Activity: Working with ultrasonic baths

DESIGNATION

Working with ultrasonic baths

DANGERS FOR HUMANS AND THE ENVIRONMENT

Dangers to humans

- Risk of hearing damage.
- Risk of explosion when exposing flammable samples to ultrasound.
- Risk of burns from the sonotrode, which can heat up to approx. 100 °C during continuous operation.

PROTECTIVE MEASURES AND RULES OF CONDUCT

Technical protective measures

- Do not remove or tamper with protective devices.
- Always use the ultrasonic processor inside the protective box.
- Only place the device on a horizontal, level and solid surface.

Organisational protective measures

- Operation only by trained persons
- Do not immerse any body parts (e.g. hands) or living beings in the bath, especially do not touch the cleaning fluid during ultrasonic cleaning.
- Do not operate the device without liquid.
- Do not fill the ultrasonic bath with hot water (more than 50°C).
- Do not put any items that contain aggressive contaminants (e.g. acids, chloride ions) into the tub.
- any flammable liquids or liquid chemicals that contain or release chloride ions (e.g. various disinfectants, household cleaners) for ultrasonic cleaning in the tub.
- Only use aggressive cleaning fluids (e.g. acids, salt solutions) in containers.
- During prolonged operation, the cleaning fluid heats up, so the temperature should be controlled when cleaning temperature-sensitive parts.
- Clean the protective box and ultrasonic bath after contamination.
- Keep the device surface and controls clean and dry.

Personal protective measures

- During prolonged continuous operation, hearing protection must be worn within a radius of 2 m.

RESPONSE TO MALFUNCTIONS

Disturbances

- In case of malfunctions, switch off the ultrasonic bath, secure it and notify the nearest supervisor.
- Never carry out repairs yourself.

WHAT TO DO IN CASE OF DISTURBANCES (continued)

Troubleshooting, repair and maintenance

- Portable electrical equipment must be inspected by a qualified expert at least once a year. The test results must be recorded in a test log.
- Repairs may only be carried out by authorized and trained personnel.
- For maintenance work, only original parts or parts that correspond to the original parts in terms of material and design may be used.
- Maintenance work may only be carried out by trained, authorized personnel (observe the manufacturer's maintenance plan).

MAINTENANCE AND REPAIR

Maintenance and repair

- Maintenance and repair work may only be carried out when the machine is at a standstill.
- Secure the device against unauthorized switching on.
- Repairs, maintenance work and inspections may only be carried out by authorized and qualified persons.

WHAT TO DO IN THE EVENT OF ACCIDENTS - FIRST AID - EMERGENCY CALL 112

- Make an emergency call.
- Initiate first aid measures.
- Trained first aiders: see first aid notice
- First aid services must be recorded in the first aid book.

CONSEQUENCES OF NON-COMPLIANCE

Injuries

- Misconduct can lead to hearing damage or burns

Property damage

- Misconduct can lead to an explosion.

Legal consequences

- Operating instructions are binding and represent written occupational safety - instructions for employees.
- Failure to comply may result in legal consequences.
- Failure to follow this instruction is a violation of the given instructions and will be punished accordingly.